

## QUICK START GUIDE (For model 2300-000)

Congratulations on your new C3 Submersible Fluorometer. We are committed to customer satisfaction. If you need assistance, technical specialists are available to answer you questions at 408-749-0994 or toll-free at 877-316-8049. This is a Quick Start Guide to help you set up the C3 Submersible Fluorometer to begin collecting data as quickly as possible.

### 1 Software Installation

1. Insert the C3 Submersible Fluorometer CD into your computer.
2. Open the C-Soft Software folder and double click on the setup icon.
3. The software will automatically be installed in the PC Program Files\C-Soft folder unless user selects alternate location.
4. There will be a prompt to restart your computer.

### 2 Connections

1. Connect the serial connector end (A) of the interface cable to the serial port on your computer.

***\*If your computer does not have a serial port use a USB to RS232 adapter.***


2. Connect the C3 Submersible Fluorometer to the 8-pin interface cable (B).

3. Connect the interface power source pin (C) to the 12 V power supply (D).

4. Plug the power supply (E) into an AC source to supply power to the C3 Submersible Fluorometer.

5. Open the C-Soft software by double clicking on the C-Soft icon.

6. If the connect icon, located in the upper corner of the C-Soft software screen, is not green, then disconnect, wait 10 seconds, and reconnect power.

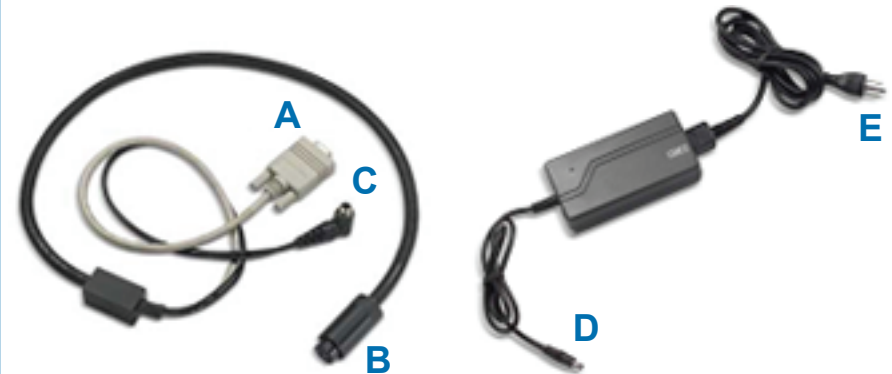

***\*The C3 Submersible Fluorometer is now ready to configure, calibrate, and initiate logging.***